

# TRACEABLE® TRIPLE-PURPOSE TIMER™ INSTRUCTIONS

## SPECIFICATIONS

Display:	3/4" high 5-digit LCD
Timing capacity:	9 hr, 59 min, 59 sec
Resolution:	1 sec
Accuracy:	0.01%
Alarm:	1 min audible/visible
Size/Weight:	3 x 2 1/2 x 1", 2 1/2 oz
Features:	Count up/stopwatch timing, countdown alarm timing, time-of-day clock, 12/24 hour formats

## CLEARING THE DISPLAY TO ZERO

Press CLEAR button to reset display to zero. In Timing mode CLEAR operates only if the timer is stopped. If the unit is timing, press START/STOP button to stop the timer before clearing display.

## CLOCK SETTING

1. Press and hold CLOCK button for 3 seconds. Bottom left cursor blinks.
2. Press AM-PM button to select the proper setting.
3. Enter correct time (all six digits) using 0-9 buttons.  
Example: If time is 4:46:18, press 0, 4, 4, 6, 1, 8. Clock is set once the last digit is entered.  
If an incorrect time is entered, press CLEAR button to clear the display.
4. Choose AM/PM or military time by pressing 12HR-24HR button located on back of unit.

## COUNT-UP TIMING

1. Press COUNT-UP button (COUNT-UP appears on display).
2. Press START/STOP button to begin count up.
3. Press START/STOP button again to stop count up (see Time-Out section).
4. With count up stopped, press CLEAR button to return the display to 0:00:00.

## COUNTDOWN TIMING

1. Press TIMER button (TIMER appears on display).
2. If necessary, CLEAR the display.
3. Enter desired countdown time using buttons 0-9.
4. Press START/STOP button to begin countdown.
5. Press START/STOP button again to stop timing (see Time-Out section).
6. When timer reaches 0:00:00, it alarms, the display flashes and begins counting up to track elapsed time since alarming.
7. When the alarm sounds, pressing the START/STOP button silences the alarm, stops the count up, and returns the display to the previously programmed time. (Must be in Timer mode to halt count up and to return display to the preset time.) **NOTE:** Alarm sounds for 1 minute then stops automatically unless turned off manually. The count up continues until turned off.

## TIME-OUT

If a pause in the countdown or count-up timing operation is required, press the start/stop button to "freeze" the display at the current reading (time-out). To resume timing at the point it was stopped, press the start/stop button again. An unlimited number of time-outs may be performed.

## ALL OPERATIONAL DIFFICULTIES

If this timer does not function properly for any reason, please replace the battery with a new high quality battery (see Battery Replacement section). Low battery power can occasionally cause any number of "apparent" operational difficulties. Replacing the battery with a new fresh battery will solve most difficulties.

## BATTERY REPLACEMENT

A faint display, incorrect display, erratic display or no display are all indicators that the battery must be replaced. On back of unit, slide open the battery cover. Remove the exhausted battery and replace with a new AAA alkaline battery. For proper placement, see drawing located inside battery compartment. Replace battery cover. Equivalent replacement battery: Control No. 1105.

## WARRANTY, SERVICE OR RECALIBRATION

For warranty, service or recalibration contact:

## CONTROL COMPANY

4455 Rex Road  
Friendswood, Texas 77546 USA  
Ph. 281-482-1714 Fax 281-482-9448  
Email sales@control3.com  
www.control3.com

Control Company is ISO 9001  
Quality-Certified by DNV  
and ISO 17025 accredited  
as a Calibration Laboratory by A2LA.

Cat. No. 5027  
Traceable® Triple-Purpose Timer Instructions  
92-5027-00  
Rev. 1 102904

Traceable® is a registered trademark of Control Company