

POST EXTENSION KIT for MULTI-TUBE VORTEXERS

1. Unplug unit from power source.
2. Remove top plate from unit.
3. Carefully place unit on its back; showing the bottom of the unit (See Fig. 1). Then remove the hex bolts from the underside of the unit. They are located under each post.
4. Turn the unit back to its upright position. Facing the side of the unit; loosen the bolts that is CLOSEST to the front of unit , 1/2 turn only (See Fig.2). This will allow the post to slide out.
5. Insert the appropriate post on each side & tighten the bolt 1/2 turn in the other direction. Make sure the groove on the longer post is facing the outside (Away) of the unit.
6. Turn unit on its back again (as in Fig. 1) re-attach the hex bolts on the underside of the unit.
7. Turn unit to its upright position.
8. Replace the top plate on unit.

Figure 1

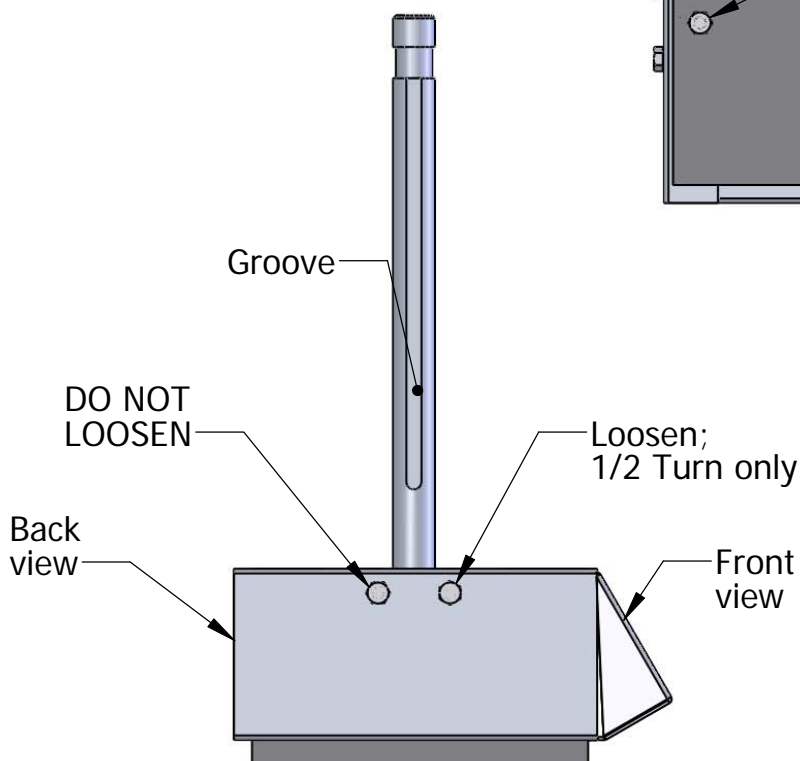
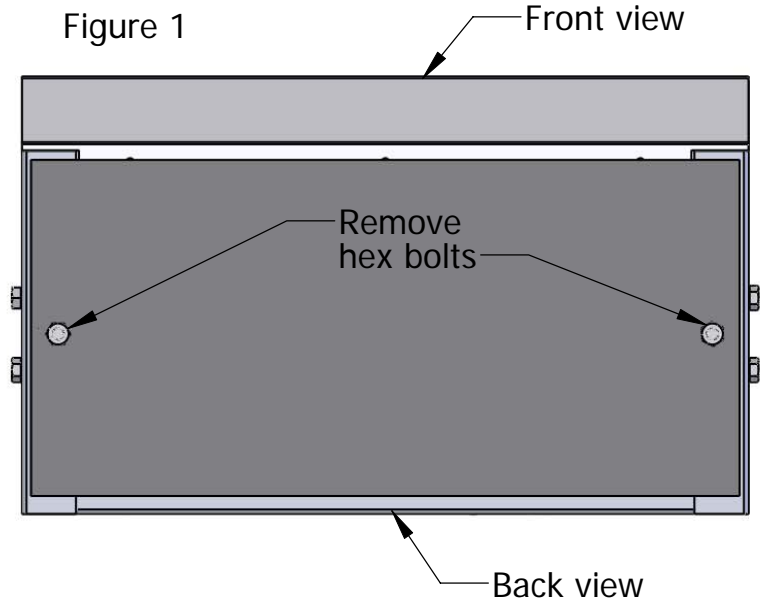


Figure 2