# HyperLadder<sup>™</sup> 50bp

Shipping: Room Temperature Batch No.: See vial Catalog numbers BIO-33054: 100 Lanes BIO-33039: 200 Lanes BIO-33040: 500 Lanes



Storage and stability: HyperLadder 50bp is shipped at room temperature. On arrival store at -20°C for optimum stability. Expiry: When stored under the recommended conditions and handled correctly, band quality is retained until the expiry date on the outer box label. Safety precautions: Please refer to the material safety data sheet for further information. Quality control specifications: HyperLadder 50bp is extensively tested quality prior to release. Notes: Research use only.

A Meridian Life Science® Company

#### Features

- Easy size determination
- 15 sharp and bright bands from 50bp 2000bp
- Ready-to-use format
- Optional mass determination

#### Description

HyperLadder 50bp (formerly HyperLadder II) is a ready-to-use molecular weight marker, especially designed for easy size determination. This ready-to-use format reduces handling steps and saves time; simply transfer HyperLadder 50bp from the vial to the gel.

HyperLadder 50bp produces a pattern of 15 regularly spaced bands, ranging from 50bp to 2000bp. To allow easy identification and orientation, the 300bp, 1000bp and 2000bp bands have the highest intensity.

Optionally the HyperLadder can be used for mass determination.

A 5x sample loading buffer is supplied for your convenience. Under no circumstances should it be used to dilute/load ladder.

### Components

Product	100 Lanes	200 Lanes	500 Lanes
HyperLadder <sup>™</sup> 50bp	500µl	2 x 500µl	5 x 500µl
5x Sample Loading Buffer	1ml	1ml	1ml


### Protocol

Vortex the tube and use  $5\mu$ l per lane.

Use 5x loading buffer for sample loading, under no circumstances should it be used for diluting the ladder. The blue dye in the loading buffer runs at approximately 4bp on a 1% gel

## **Associated Products**

Product	Pack size	Cat. No.
Agarose	100g	BIO-41026
Agarose Tablets	150g	BIO-41028

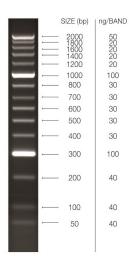

#### Notes

- 1. This product insert is a declaration of analysis at the time of manufacture
- 2. HyperLadder is a trademark of Bioline
- 3. No licence is conveyed in respect of US 6,680,378
- 4. Research Use Only

Bioline Reagents Ltd UNITED KINGDOM

Tel: +44(0)20 8830 5300 Fax: +44 (0)20 8452 2822

PI-50095 V3

Fax: +1 508 880 8993

Bioline USA Inc.

Tel: +1 508 880 8990

USA

Bioline GmbH GERMANY

Tel: +49(0)33 7168 1229 Fax: +49 (0)337168 1244 Bioline (Aust) Pty. Ltd AUSTRALIA

Tel: +61 (0)2 9209 4180 Fax: +61 (0)2 9209 4763