

ProSieve™ Color Protein Markers

Volume: 500 μ l Cat. No. 50550
Volume: 100 μ l Cat. No. 50552

Store: -20°C

For best results, lyophilized material should be stored at -20°C and used within five years from date of manufacture (DOM).

Stable for three months once reconstituted. Do not freeze/thaw reconstituted vials of ProSieve™ Color Marker more than five times.

Recommended procedure

First Time Use

1. Add 500 μ l water (Cat. No. 50550) or 100 μ l water (Cat. No. 50552).
2. Warm briefly to 37°C and mix well.
3. Load 10 μ l per well on standard polyacrylamide mini-gels. ProSieve™ Color Protein Markers are ideal for use with PAGEr™ Gold Precast Gels.
4. Aliquot and store at -20°C.

Subsequent Use

Warming aliquots to 37°C is not required for subsequent use of reconstituted ProSieve™ Color Protein Marker.

For Research Use Only.

All trademarks herein are marks of the Lonza Group or its affiliates.

© 2007 Lonza Rockland, Inc.
All rights reserved.