



## Quick-ITS<sup>™</sup> Plus NGS Library Prep Kit

Fastest, normalization-free ITS rRNA library prep

### **Highlights**

- The most streamlined NGS kit with only 30 minutes of hands-on time • for 96 samples.
- 100% automation ready with only a single PCR step and without the • need for normalization.
- Real-time PCR enables absolute microbial number . copy quantification.

Catalog Numbers: D6424, D6426



Scan with your smart-phone camera to view the online protocol/video.







## **Table of Contents**

Product Contents	01
Specifications	
Product Description	03-04
Protocol	05-09
Before Starting	
(I) 1-Step PCR	05-07
(II) Pooling by Equal Volume	07
(III) Final Library Clean-up	
(IV) Library Quantification	
Appendices	10-12
Absolute Quan tification of Total ITS Copy Number by Real-Time PCR Improving Normalization by Fine-Tuning	10
Pooling Volumes based on qPCR	11
Index Primer Sets	11
Removal of PCR Inhibitors from DNA Composition of ZymoBIOMICS <sup>™</sup>	11
Microbial Community DNA Standard	12
Illumina MiSeq <sup>®</sup> Sample Sheet Setup	12
Ordering Information	13
Complete Your Workflow	14
Troubleshooting Guide	15
Notes	16
Guarantee	17

### **Product Contents**

<i>Quick</i> -ITS <sup>™</sup> Plus NGS Library Prep Kit	D6424 (96 rxns.)	D6426 (24 rxns.)	Storage Temp.
Equalase <sup>™</sup> qPCR Premix	1 ml	400 µl	-20°C
ZymoBIOMICS <sup>™</sup> DNase/RNase Free Water	1 ml	1 ml	Room Temp.
ZymoBIOMICS <sup>™</sup> Microbial Community DNA Standard (50 ng)	10 µl	10 µl	-20°C
ZymoBIOMICS <sup>™</sup> 16S/ITS qPCR Standard	10 µl	10 µl	-20°C
ITS Index Primer Set <sup>1</sup>	A, 10 μl each (plate)	Z, 20 µl each (tube)	-20°C
PCR Inactivation Solution	100 µl	100 µl	Room Temp.
Select-a-Size MagBead Concentrate <sup>2</sup>	30 µl	30 µl	4-8°C
Select-a-Size MagBead Buffer <sup>2</sup>	1 ml	1 ml	4-8°C
DNA Wash Buffer	6 ml	6 ml	Room Temp.
Magnetic Rod	4	4	-
Instruction Manual	1 pc	1 pc	-

<sup>1</sup> ITS Index Primer Set A contains the Index Primers ITSR ZT725-ZT736 and ITSF ZT517-ZT524 already combined together to reach 10 µl total (5 µl of i7 index and 5 µl of i5 index) in a 96-well microplate. ITS Index Primer Set Z contains the Index Primers ITSR ZT725-ZT730 and ITSF ZT517-ZT520 in tubes. The barcodes of each index primer are distinct from one another by at least 5 bp to boost the accuracy of demultiplexing. Note that the i5 bases remain the same on the sample sheet, while thei7 bases are entered as the reverse complement on the sample sheet. See Appendix F for additional support and a template for the Illumina MiSeq<sup>®</sup> sample sheet.

<sup>2</sup> The Select-a-Size MagBead Concentrate and Buffer are shipped at room temperature but should be stored at 4-8°C upon receipt.

### **Specifications**

- **Sample Input –** Purified microbial DNA (≤100 ng), free of PCR inhibitors<sup>1</sup>.
- ITS Primer Sequences (adapters not included)<sup>2</sup> ITS3f (GCATCGATGAAGAACGCAGC, 20 bp), ITS4r (TCCTCCGCTTATTGATATGC, 20 bp).
- Index Primers Dual index (barcodes) to uniquely label samples<sup>3</sup>.
- Barcode Sequences <u>10 bp</u> barcodes are available for download by visiting the Documentation section of the D6424/D6426 Product Page at <u>www.zymoresearch.com</u>.
- **Amplicon Size** The final amplicon size after 1-Step PCR (targeted amplification and barcode addition) is ~480 bp.
- Sequencing Platform Illumina MiSeq<sup>®</sup> without the need to add custom sequencing primers. We recommend the MiSeq<sup>®</sup> Reagent Kit v3 (600-cycle). For assistance with sample sheet setup, see Appendix F.
- Equipment Needed (user provided) Microcentrifuge, plate spinner (centrifuge), 96-well real-time quantitative PCR system (SYBR Green compatible, recommended) or standard PCR system, and 96-well real-time PCR plates.

<sup>1</sup> DNA that contains potent PCR inhibitors such as polyphenolics, humic/fulvic acids, tamins, melanin, etc. can be quickly cleaned using the OneStep<sup>™</sup> PCR Inhibitor Removal Kit. See Appendix D for additional information. 2 Primers from White, T.J. et al. (1990). Amplification and Direct Sequencing of Fungal Ribosomal RNA Genes for Phylogenetics. PCR Protocols and Applications, pp.315-322.

<sup>3</sup> For projects with more than 96 samples, please contact Zymo Research at <u>oemorders@zymoresearch.com</u> for additional indexing solutions.

### **Product Description**

ITS rRNA gene sequencing is a routine technique for mycobiome (fungal) composition profiling. Compared to shotgun metagenomics sequencing, ITS rRNA gene sequencing is more cost-effective and more robust; it generally requires less input DNA and is less impacted by the presence of non-microbial DNA. However, ITS rRNA gene sequencing has its own challenges. One major challenge is the formation of PCR chimeric sequences, which are artificial sequences resulting from the recombination of two or more PCR templates. Moreover, common ITS library preparation protocols have not been optimized to be cost-effective for large-scale applications.

The **Quick-ITS**<sup>™</sup> **Plus NGS Library Prep Kit** is the fastest and simplest library prep method for high-throughput ITS rRNA sequencing. Distinguishing features of the kit are described below.

Fastest ITS rRNA Workflow. The Quick-ITS<sup>™</sup> Plus NGS Library Prep Kit utilizes a single qPCR/PCR for combined targeted amplification and barcode addition using specially designed primers. After pooling by equal volume, a single clean-up of the final library is performed, rather than multiwell magnetic bead clean-ups. Additional library quantification analysis such as TapeStation<sup>®</sup> analysis or gel electrophoresis are not necessary. With this workflow, the hands-on time of ITS library preparation is reduced to only 30 minutes (Figure 1).

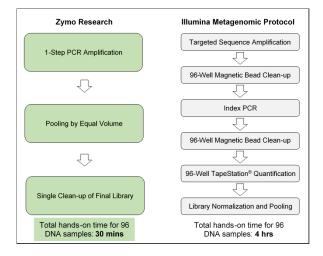


Figure 1. Quick-ITS<sup>™</sup> Plus NGS Library Prep Kit workflow versus the Illumina Metagenomic Protocol. Total hands-on time calculations are based on the preparation of 96 DNA samples.

**Normalization Free.** There is no need to control sample input or PCR cycles. The workflow auto-normalizes and produces similar amounts of sequencing reads across all samples regardless of different DNA inputs (Figure 2). Just pool by equal volumes!

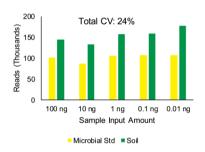


Figure 2. The Quick-ITS<sup>™</sup> Plus NGS Library Prep Kit results in similar amounts of reads across different input amounts without normalization. 0.01-100 ng ZymoBIOMICS<sup>™</sup> Microbial Community Standard and soil DNA were used as inputs. Libraries were pooled by equal volumes (2 µl each) without further normalization and sequenced using the MiSeq<sup>®</sup> Reagent Kit v3 (600-cycle). The CV (coefficient of variation) is the ratio of the standard deviation to the mean with lower values corresponding to less dispersion around the mean.

**High Quality NGS Library.** The workflow has been optimized to minimize amplification bias (Figure 3). PCR chimera formation is maintained below 10% without the need to PCR cycles (Figure 4).

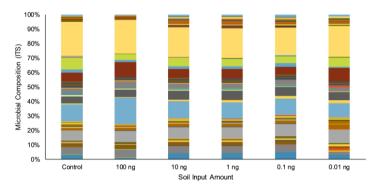


Figure 3. Compatible with wide sample input range. Fungal composition profiles (species level shown) are accurate from 100 ng down to 0.01 ng using soil DNA input and agree with the control (Illumina).

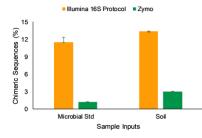


Figure 4. Up to 70% chimera reduction without controlling PCR cycles. Equal amounts of ZymoBIOMICS<sup>™</sup> Microbial Community Standard and soil DNA were used as inputs and amplified for 42 cycles using 1-step PCR and Equalase<sup>™</sup> compared to using the Illumina Metagenomic Protocol. Chimeric sequences were predicted with UCHIME (https://www.drive5.com/uchime).

### Protocol

#### **Before Starting**

- ✓ Sample Quantity Requirement. To ensure color balance in index sequencing, a minimum of 9 samples per run is recommended using the index primers (ZT517-ZT519, ZT725-ZT727).
- Input DNA Guidelines. All DNA samples should be free of PCR inhibitors<sup>1</sup>.

#### Section 1: 1-Step PCR

1. Set up a master mix<sup>2</sup> according to the table below:

Component	Volume/Reaction
Equalase <sup>™</sup> qPCR Premix	10 µl
ZymoBIOMICS <sup>™</sup> DNase/RNase Free Water	4 µl
Total	14 µl

- For each reaction, add 14 µl of the master mix to the appropriate wells of a 96-well real-time PCR plate. A sample of the plate setup can be found on the next page on the Plate Setup Guide<sup>3</sup>.
- From ITS Index Primer Set<sup>4</sup>, pierce the foil and add 4 μl of the appropriate Index Primer ITR ZT7XX and Index Primer ITSF ZT5XX combination to the proper wells of the PCR plate as indicated in the diagram on the next page:

(Continued on next page.)

<sup>1</sup> DNA that contains potent PCR inhibitors such as polyphenolics, humic/fulvic acids, tarnins, melanin, etc. can be quickly cleaned using the OneStep<sup>™</sup> PCR Inhibitor Removal Kit. See Appendix D for additional information. 2 Prepare a master mix to include both a positive and negative control.

<sup>3</sup> The Plate Setup Guide is available for download by visiting the Documentation section of the D6424/D6426 Product Page at www.zymoresearch.com.

<sup>4</sup> If using ITS Index Primer Set Z, add 2 µl of i7 index primer and 2 µl of i5 index primer from the appropriate tubes.

Index Primers ITSR ZT7xx														
			ZT725	ZT726	ZT727	ZT728	ZT729	ZT730	ZT731	ZT732	ZT733	ZT734	ZT735	ZT736
			1	2	3	4	5	6	7	8	9	10	11	12
	ZT517	А	S1	S9	\$17	S25	\$33	S41	S49	\$57	S65	\$73	S81	S89*
×	ZT518	в	S2	S10	S18	S26	\$34	S42	\$50	S58	S66	S74	S82	\$90*
ZT5xx	ZT519	с	S3	S11	S19	\$27	\$35	S43	S51	S59	S67	S75	S83	S91*
TSF	ZT520	D	S4	\$12	S20	S28	\$36	S44	\$52	\$60	S68	S76	S84	S92*
ers l	ZT521	E	S5	S13	S21	S29	\$37	S45	S53	S61	S69	S77	S85	S93*
Index Primers ITSF	ZT522	F	S6	S14	\$22	\$30	S38	S46	S54	S62	\$70	S78	S86	\$94*
lex F	ZT523	G	S7	\$15	S23	\$31	\$39	S47	\$55	S63	S71	S79	S87	POS**
Ē	ZT524	н	S8	S16	S24	S32	S40	S48	S56	S64	S72	S80	S88	NEG***
<ul> <li>*** POS: The ZymoBIOMICS<sup>™</sup> Microbial Community DNA Standard<sup>1</sup> (included in kit) as a positive control.</li> <li>*** NEG: A no template control as a negative control.</li> <li>4. (Optional): If absolute quantification by real-time PCR is desired, add 2 µl of the serially diluted qPCR standard to the 6 wells highlighted above; S89-S94. Refer to Appendix A for more details.</li> <li>5. Add 2 µl of your DNA samples to individual wells. Include a positive and negative control in the plate.</li> <li>6. Apply an adhesive PCR plate seal. Mix the plate on a plate shaker and centrifuge in a plate spinner<sup>2</sup>.</li> <li>7. Place plate in a real-time thermocycler<sup>3</sup> and run the program shown</li> </ul>														
TemperatureTime $95^{\circ}C$ 10 min $95^{\circ}C$ 30 sec $55^{\circ}C$ 30 sec $72^{\circ}C$ 3 minPlate read- $4^{\circ}C$ Hold(Continued on next page.)														

<sup>1</sup> The composition of the microbial standard can be found in Appendix E.

<sup>2</sup> PCR reactions can be pipette mixed if a plate shaker is not available.

<sup>3</sup> A real-time thermocycler is recommended as it enables QC of the library prep of all wells and absolute quantification as shown in Appendix A. A non-quantitative system can be used if absolute quantification is not needed.

- Once the samples have cooled to 4°C, stop the program. Centrifuge plate in a plate spinner to collect condensation in wells and place plate on ice. Proceed to <u>Section 2</u>, or store plate at ≤-20°C for later use.
- 9. Monitor and QC the library preparation when running the reaction on a real-time thermocycler<sup>1</sup>.
  - a. For example, a sample that is expected to amplify and shows little or no amplification may indicate an error in the reaction setup (See the Troubleshooting Guide).
  - b. The negative control should not amplify before 35 cycles<sup>2</sup>. Earlier amplification of negative control might indicate process contaminations.
  - c. An example of qPCR amplification with controls is shown in Figure 5 below.

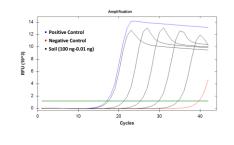


Figure 5. qPCR Amplification Example with Positive and Negative Controls. Serial dilution of soil DNA (black) from 100 to 0.01 ng to were amplified on a Bio-Rad CFX96<sup>™</sup> Real-Time PCR Detection System. The positive (blue) amplified at 17.01 and negative (red) amplified at 39.60. Baseline threshold was set at 1200 RFU.

#### Section 2: Pooling by Equal Volume

Add 50 µl of **PCR Inactivation Solution** into a new microcentrifuge tube. Pool equal volumes (5 µl) of PCR products from each well of the plate from <u>Section 1</u> into the tube<sup>3</sup> and mix well. Skip the wells of S89-S94 if they are used for qPCR standards. Proceed to <u>Section 3</u>.

<sup>1</sup> If real-time PCR was not used, after amplification perform PCR cleanup for a few samples plus positive and control. Analyze on a TapeStation<sup>®</sup> to confirm correct amplicon size (~480 bp).

<sup>2</sup> The PCR program runs for 42 cycles, so it is normal to see some amplification from the negative control. The negative control should be sequenced together with other samples. If appropriate for your project, the taxa from the negative control can be subtracted from the analysis.

<sup>3</sup> There is no need for additional normalization procedures. To obtain better normalization and adjust pooling volumes, See Appendix B.

#### Section 3: Final Library Clean-up

- Equilibrate the Select-a-Size MagBead Buffer to room temperature (15-30°C). Add 30 µl of Select-a-Size MagBead Concentrate to the 1 ml Select-a-Size MagBead Buffer. Resuspend the magnetic particles by vigorously shaking until homogenous.
- Add Select-a-Size MagBeads to the pooled library from <u>Section</u> <u>2</u> at a ratio of 0.8x volume. For example, add 400 µl of Select-a- Size MagBeads to 500 µl of the pooled library and PCR Inactivation Solution mixture.
- 3. Mix thoroughly by pipetting or vortexing until homogenous. Incubate for 5 minutes at room temperature.
- 4. Place the sample on a magnetic rack<sup>1</sup> and incubate for 3-10 minutes at room temperature, or until the magnetic beads have fully separated from solution.
- 5. Once the beads have cleared from solution, remove and discard the supernatant<sup>2</sup>.
- While the beads are still on the magnetic rack, add 1 ml of DNA Wash Buffer. Remove and discard the supernatant. <u>Repeat this</u> <u>step</u>.
- 7. While the beads are still on the magnetic rack, aspirate out any residual buffer with a 10 µl pipette tip.
- 8. Remove tube from the magnetic rack and keep the cap open for 3 minutes at room temperature to dry the beads.
- Add 10-100 µl<sup>3</sup> of ZymoBIOMICS<sup>™</sup> DNase/RNase Free Water to the beads and pipette mix thoroughly. Incubate at room temperature for 2 minutes.
- 10. Place the sample on a magnetic rack and incubate for 1 minute at room temperature, or until the magnetic beads have fully separated from eluate.
- 11. Transfer supernatant to a clean microcentrifuge tube. Proceed to <u>Section 4</u>.

<sup>1</sup> Alternatively, the provided Magnetic Rod can be used.

<sup>2</sup> Avoid aspirating any beads when removing the supernatant. To best prevent this, leave 2-5 µl of liquid behind. 3 If pooling fewer than 10 samples, use 10 µl for elution.

### Section 4: Library Quantification

Use a fluorescence-based method (Qubit<sup>®</sup> dsDNA HS Assay Kit recommended) to quantify the final library. Using a final amplicon size of 480 bp, convert ng/µl to nM using the equation below.

 $\frac{\text{concentration in } ng/ul}{660 \ g/mol \ x \ average \ library \ size \ in \ bp} \ x \ 10^6 = \text{concentration in } nM$ 

For example: 20 ng/µl DNA of the final library is equivalent to 63.13 nM.

If preferred, a qPCR-based method for quantification may be used<sup>1</sup>.

#### **DNA Fragment Analysis (Not Required)**

If a fragment analyzer (e.g. TapeStation<sup>®</sup>) is used to analyze the final library, there may be a lack of a tight band at ~480 bp. Because of the library prep design, some library products have run through additional PCR cycles and might not anneal well. They are perfectly fine for sequencing; which already denatures double-stranded DNA into single strands.

#### This is your final ITS library.

The ultra-pure pooled library DNA is now ready for use or storage at ≤-20°C. Refer to platform-specific guidelines for preparation for sequencing. <u>No custom sequencing primers are needed</u>.

### Illumina MiSeq<sup>®</sup> Setup:

The MiSeq<sup>®</sup> Reagent Kit v3 (600-cycle) with 15% PhiX spike-in is recommended. See Appendix F for assistance with sample sheet setup. <u>Remember to set the index size to 10 bp.</u>

1 For example, KAPA® Library Quantification Kit or similar.

## Appendices

# Appendix A: Absolute Quantification of Total ITS Copy Number by Real-Time PCR

Absolute quantification is determined with a real-time PCR standard curve using serially diluted amounts of the ZymoBIOMICS<sup>™</sup> 16S/ITS qPCR Standard.

- Dilute the ZymoBIOMICS<sup>™</sup> 16S/ITS qPCR Standard (7.5x10<sup>6</sup> copies/µl) to 7.5x10<sup>6</sup> copies/µl, 7.5x10<sup>4</sup> copies/µl and 7.5x10<sup>2</sup> copies/µl.
- Add 2 µl of the standard and its dilutions to six wells (each in duplicate in positions S89-S94; see plate layout on Page 6) on the same plate that contains unknown samples for library prep.
- After 1-Step PCR, plot the concentration (ITS copies/µI) of the standards against the Ct values to produce a standard curve. See Figure 6 and 7 for an example.
- 4. Calculate the total ITS copy number of unknown samples based on their Ct value with the standard curve accordingly<sup>1</sup>.

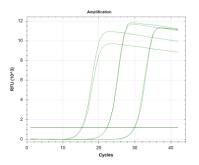


Figure 6. qPCR of Diluted Amounts of 165/ITS Standard. The standard was amplified on a Bio-Rad CFX96<sup>™</sup> Real-Time PCR Detection System. Baseline threshold was set at 1200 RFU.

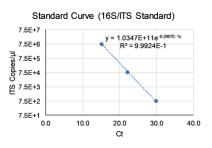


Figure 7. Absolute Quantification using a Standard Curve. Ct can be correlated to the total ITS copies based on the regression.

<sup>1</sup> Excel template for the Absolute Quantification of Total ITS Copy Number is available for download by visiting the Documentation section of the D6424/D6426 Product Page at <a href="http://www.zymoresearch.com">www.zymoresearch.com</a>.

### Appendix B: Improving Read Distribution by Fine-Tuning Pooling Volumes based on qPCR

Typically, pooling samples by equal volumes will result in sufficient normalization and similar number of reads per sample. To obtain even better normalization, the pooling volumes can be adjusted based on Ct values obtained.

First, adjust the fluorescence threshold/cutoff to a certain value so that the positive control<sup>1</sup> has a Ct of 17. Use the following equation to calculate the pooling volume of each well based on its Ct: <u>Volume (µl) =  $0.00104x^2 - 0.0965x + 3.736$ ; x = Ct</u>. For example, the pooling volumes for Ct values of 15, 20, and 30 are 2.52 µl, 2.22 µl, and 1.78 µl respectively. This strategy can further reduce the CV (coefficient of variation) of reads assigned to samples by 3-6%.

### Appendix C: Index Primer Sets

To accommodate sequencing projects of various sizes, Zymo Research offers primer sets that can uniquely barcode more than 96 samples. The barcodes of each index primer are distinct from one another by at least 5 bp to boost the accuracy of demultiplexing. For projects that require indexes for >96 samples, please contact Zymo Research at <u>oemorders@zymoresearch.com</u> for a custom solution.

### Appendix D: Removal of PCR Inhibitors from DNA

The input DNA samples for the Quick-ITS<sup>™</sup> Plus NGS Library Prep Kit must be free of PCR inhibitors such as polyphenolics, humic/fulvic acids, tannins, melanin, etc. To further remove PCR inhibitors from purified DNA samples, Zymo Research recommends performing a one-step cleanup with the OneStep<sup>™</sup> PCR Inhibitor Removal Kit. Additional information can be found by visiting the D6030 Product Page at <u>www.zymoresearch.com</u>.

<sup>1</sup> The positive control is the sample that uses 2  $\mu l$  of the ZymoBIOMICS  $^{\circ}$  Microbial Community DNA Standard.

### <u>Appendix E: Composition of ZymoBIOMICS<sup>™</sup> Microbial Community</u> <u>DNA Standard (50 ng)</u>

The **ZymoBIOMICS<sup>™</sup> Microbial Community DNA Standard (50 ng)** is a mixture of genomic DNA extracted from pure cultures of eight bacterial and two fungal strains. Genomic DNA from each culture is quantified before mixing. The ZymoBIOMICS<sup>™</sup> Microbial Community DNA Standard allows for assessment of bias from library preparation, sequencing, and bioinformatics analysis. More information about the standard can be found below. For more detail, see D6305 Instruction Manual.

	Theoretical Composition (%)					
Species	Genomic DNA	16S Only	16S & 18S	Genome Copy	Cell Number	
Pseudomonas aeruginosa	12	4.2	3.6	6.1	6.1	
Escherichia coli	12	10.1	8.9	8.5	8.5	
Salmonella enterica	12	10.4	9.1	8.7	8.8	
Lactobacillus fermentum	12	18.4	16.1	21.6	21.9	
Enterococcus faecalis	12	9.9	8.7	14.6	14.6	
Staphylococcus aureus	12	15.5	13.6	15.2	15.3	
Listeria monocytogenes	12	14.1	12.4	13.9	13.9	
Bacillus subtilis	12	17.4	15.3	10.3	10.3	
Saccharomyces cerevisiae	2	NA	9.3	0.57	0.29	
Cryptococcus neoformans	2	NA	3.3	0.37	0.18	

#### Figure 8: Microbial Composition

### Appendix F: Illumina MiSeq<sup>®</sup> Sample Sheet Setup

A template for the Illumina MiSeq<sup>®</sup> sample sheet is available for download by visiting the Documentation section of the Product Page at <u>www.zymoresearch.com</u>. Fill in the project and sample information in the highlighted fields, then save the file in comma-separated values (CSV) format for use with the Illumina MiSeq<sup>®</sup>.

# **Ordering Information**

Product Description	Catalog No.	Size
<i>Quick</i> -ITS <sup>™</sup> Plus NGS Library Prep Kit	D6424 D6426	96 rxns. 24 rxns.

Individual Kit Components	Catalog No.	Amount
ZymoBIOMICS <sup>™</sup> DNase/RNase Free Water	D4302-5-10	10 ml
ZymoBIOMICS <sup>™</sup> Microbial Community <u>DNA</u> Standard (200 ng)	D6305	200 ng
ZymoBIOMICS <sup>™</sup> Microbial Community <u>DNA</u> Standard (2000 ng)	D6306	2000 ng

## **Explore Other Microbiome Products**

- DNA/RNA Shield<sup>™</sup> and Collection Devices

   1X Reagent #R1100
   For sample lysis and stabilization of DNA/RNA

   2X Concentrate #R1200
   Reagent concentrate (2X) for use with liquids at 1:1 ratio

   Fecal Collection Tube #R1101
   15 mL container (prefilled with 9 mL DNA/RNA Shield<sup>™</sup>). Direct collection of up to 1g or 1 mL stool

   Collection Tube w/ Swab #R1106
   12 x 80 mm screwcap container filled with 1 mL DNA/RNA Shield<sup>™</sup> and sterile swab for specimen collection
- To collect and transport samples at ambient temperatures:

✓ Unbiased and inhibitor-free DNA and RNA extraction (high-throughput and automatable) for microbial profiling:

	ZymoBIOMICS <sup>™</sup> DNA and RNA Kits	
6	DNA Miniprep #D4300	Up to 25 µg DNA
	DNA Microprep #D4301	Up to 5 µg DNA
	MagBead DNA #D4302	Automatable (Tecan, Hamilton, Kingfisher, etc.)
	96-Well DNA #D4309	Spin-plate
	DNA/RNA Miniprep Kit #R2002	Up to 100 μg DNA/RNA

 Microbial standards and references for profiling quality control, ben chmarking, positive controls, and to assess performance of entire microbiomic/metagenomic workflows:

	ZymoBIOMICS <sup>®</sup> Standards and Refer	ence Materials
	Microbial Community Standard #D6300	Contains 8 bacteria and 2 yeasts for QC and method optimization
	Microbial Community DNA Standard #D6305	Contains 8 bacteria and 2 yeasts DNA for bioinformatics optimization
	Gut Microbiome Standard #D6331	Contains 21 different human gut strains for method benchmarking
	Fecal Reference with TruMatrix <sup>™</sup> Technology #D6323	Contains real human fecal material for benchmarking and improved data reproducibility

# **Troubleshooting Guide**

Problem	Possible Causes and Suggested Solutions
Background Contamination	Workspace contamination: - Clean workspace, microcentrifuge, and pipettes with 10% bleach routinely to avoid contamination.
	<ul> <li>Use of kit in exposed environment without proper filtration can lead to back ground contamination. Check pipettes, pipette tips, microcentrifuge tubes, workspace, etc. for contamination</li> </ul>
	<ul> <li>Make sure all reagent tubes and bottles are properly sealed for storage.</li> <li>Use of these outside a clean room or hood can result in contamination.</li> </ul>
Loss of Volume during PCR	Adhesive seal: - A loosened adhesive seal on the PCR plate can lead to sample evaporation. Ensure that the plate seal is secure on every well during targeted sequence amplification. Lid pressure:
	<ul> <li>Inconsistent lid pressure. Ensure that the lid pressure on the real-time quantitative PCR instrument is consistent over the PCR plate according to the manufacturer's recommendation.</li> </ul>
Unexpected or No	Sample with high microbial DNA concentration:
Amplification of DNA Sample During PCR Program in <u>Section 1</u>	<ul> <li>Reaction setup error. A sample that is expected to amplify but shows little or no amplification during the PCR program in <u>Section 1</u> may indicate an error in the reaction setup. Use a new aliquot of the sample and repeat <u>Section 1</u>.</li> </ul>
	Sample with low microbial DNA concentration:
	<ul> <li>Check negative control. A sample with little microbial DNA may not amplify before the negative control. Either use more concentrated DNA or use more DNA volume during reaction setup.</li> </ul>
	Abnormal qPCR curves:
	<ul> <li>Proceed as normal. Abnormal qPCR amplification curves may occur and this is normal performance. This is usually a slight dip in RFU (forming a small "hump").</li> </ul>
Diminished Amplicon Bands in Library Analysis	No single amplicon peak and/or high background: - Proceed as normal. There may be a lack of a single band and/or high background if using TapeStation <sup>®</sup> or similar methods to determine amplicon size. This is normal and is part of the library prep design. Do not use this sizing and quantification data. To properly quantify library, use a fluorescence-based method and calculation in <u>Section 4</u> .

For technical assistance, please contact 1-888-882-9682 or email tech@zymoresearch.com

### Notes




### 100% satisfaction guarantee on all Zymo Research products, or your money back.

Zymo Research is committed to simplifying your research with quality products and services. If you are dissatisfied with this product for any reason, please call 1(888) 882-9682.

Integrity of kit components is guaranteed for up to one year from date of purchase. Reagents are routinely tested on a lot-to-lot basis to ensure they provide the highest performance and reliability.

This product is for research use only and should only be used by trained professionals. It is not for use in diagnostic procedures. Some reagents included with this kit are irritants. Wear protective gloves and eye protection. Follow the safety guidelines and rules enacted by your research institution or facility.

<sup>™</sup> Trademarks of Zymo Research Corporation, <sup>®</sup> Registered Trademarks of Zymo Research Corporation Other trademarks: Bio-Rad CFX96<sup>™</sup> Real-Time PCR Detection System (Bio-Rad Laboratories, Inc.), MiSeq<sup>®</sup> (Illumina, Inc.), TapeStation<sup>®</sup> (Agilent Technologies, Inc.), Qubit<sup>®</sup> (Thermo Fisher Scientific). Kapa<sup>®</sup> (Roche), TruMatrix<sup>™</sup> (The BioCollective).



### The **BEAUTY** of **SCIENCE** is to Make Things **SIMPLE**<sup>®</sup>





