

Mini-Alarm-Timer/ Stopwatch Instructions

SPECIFICATIONS

- Display:** 1/2"-High LCD showing minutes and seconds
- Accuracy:** 0.01%
- Timing**
- Capacity:** Count down – 1 second to 99 minutes, 59 seconds Count-up – 1 second to 59 minutes, 59 seconds.
- Alarm:** Audible alarm sounds for one minute
- Functions:** Countdown timing and count-up (stopwatch) timing
- Attachments:** Clip, magnetic back, and stand
- Size:** 2 1/2" x 2 1/4" x 1/2"
- Weight:** 1.5 oz

COUNTDOWN TIMING

1. Press the STOP/RESET button to set timer to zero. The display should read 00:00.
2. Press the numbers keys 0 through 9 to set desired minutes and seconds. (Example: to set 12 minutes 38 seconds, press 1, 2, 3, 8.)

Note: To time in seconds only, press just two of the number keys. (Example: To set 15 seconds, press keys 1 and 5.)

3. If the wrong time has been entered, press the STOP/RESET button to return the display to 00.00.
4. Once the desired time is on the display, press the START button to begin countdown timing.
5. When the timer reaches 00:00, an alarm will sound. The alarm will sound for 60 seconds and then turn off automatically. To turn the alarm off manually, press any button or number key.

Note: The display will read 00.00 during the 60 second alarm sounding time and after the alarm is turned off.

6. During the countdown process, the timer may be paused by pressing the STOP/RESET button one time and restarted by pressing the START button.

Note: Pressing the STOP/RESET button twice, will clear the display and return it to 00.00.

STOPWATCH (COUNT-UP) TIMING

1. Press the STOP/RESET button to set the timer to zero. The display should read 00.00.
2. Press the START button once to begin count-up timing.

Note: Timer will count up to 59 minutes, 59 seconds and then automatically reset itself to zero – the display will read 00.00 for one second and then start counting up

3. During the count-up process, the timer may be paused by pressing the STOP/RESET button one time and restarted by pressing the START button.

Note: Pressing the STOP/RESET button twice, will clear the display and return it to 00.00.

4. To end the count-up process at any time, press the STOP/RESET button twice.

TIMEOUT

The timing process may be stopped at any time during count-up or countdown. Press the STOP/RESET button one time. Timing may be resumed by pressing the START button.

Note: Pressing the STOP/RESET button twice, will clear the display and return it to 00.00.

All Operational Difficulties

If this timer does not function properly for any reason, please replace the battery with a new high quality battery (see "Battery replacement" section). Low battery power can occasionally cause any number of "apparent" operational difficulties. Replacing the battery with a new fresh battery will solve most difficulties.

Battery Replacement

An incorrect display, no display or operational difficulties indicate that the battery should be replaced. Use a coin to open the battery cover on the back of the timer (turn the cover approximately 1/8 of a turn counter-clockwise). Equivalent battery replacements are: RAYOVAC RW42, DURACELL D357, GPA76, and EVEREADY 357. Insert the battery with the positive side facing you. Replace the battery cover. Replacement battery Cat. No. 1039.

WARRANTY, SERVICE, RECALIBRATION

For warranty, service or recalibration contact:

CONTROL COMPANY

4455 Rex Road
Friendswood, Texas 77546 USA
Ph. 281-482-1714 Fax 281-482-9448
E-mail sales@control3.com
www.control3.com

Control Company is ISO 9001 Quality-Certified by DNV and ISO 17025 accredited as a Calibration Laboratory by A2LA.

Cat. No. 5020

Traceable® Mini-Alarm-Timer/Stopwatch Instructions
92-5020-00
Rev. 0 040705

Traceable® is a registered trademark of Control Company

© 2005 Control Company