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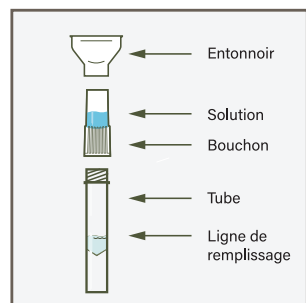
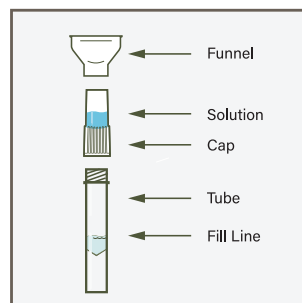
Fill the tube with saliva to the black wavy line.

Fill the tube until your saliva (not including bubbles) is at or just above the wavy line. **DO NOT OVERFILL.**

Replace the funnel with the fluid cap.

Remove the funnel from the tube. Screw on the enclosed cap **TIGHTLY** to release the solution that will stabilize the DNA in your saliva.

Firmly screw cap down to release solution and seal tube.

You will know it works when the blue solution from the cap is released into the tube. Firmly tighten cap to assure the cap and tube is completely sealed.

Shake the tube for at least five seconds.

This will ensure your sample mixes thoroughly with the stabilizing solution.

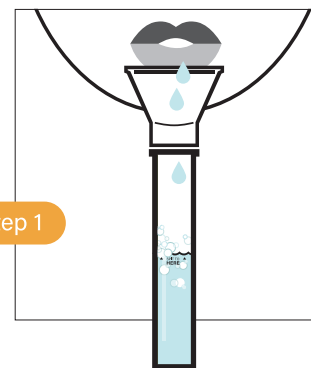
Cap choking hazard. Keep out of reach of children. Wash with water if liquid solution comes in contact with eyes or skin. **DO NOT INGEST LIQUID.**

This kit is designed for the collection of human saliva samples.


Made in USA, Patent Pending

For research use only. Not for use in diagnostic procedures.


SPECTRUMDNA™
A Division of SPECTRUM SOLUTIONS, LLC

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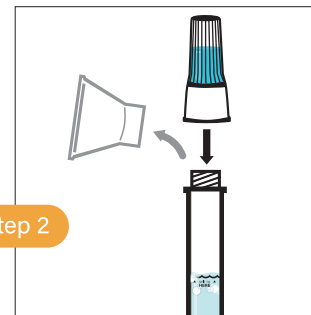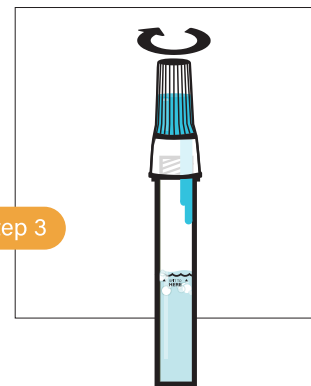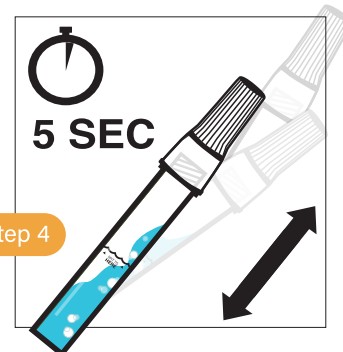
Label Legend:
I Consult instructions for use
U Use by date (Collect saliva by)
IVD In vitro diagnostic medical device
REF Catalog number
CE CE Marking
⚠ Caution, consult instructions for use
⬇ Temperature limit
IC/IMP Authorized representative
■ Manufacturer
LOT Batch code