


TRACEABLE® FOOD/WATERPROOF/PIERCING/DROP-PROOF THERMOMETER INSTRUCTIONS

SPECIFICATIONS:

Display: 1/4" LCD
Range: -58 to 536 °F / -50 to 280 °C
Resolution: 0.1° (-20 to 200), 1° outside of range
Accuracy: **Cat. No. 4369** - $\pm 0.4^{\circ}\text{C}$ at tested points
Cat. No. 4370 - $\pm 1.0^{\circ}\text{C}$
Sampling Rate: 1 second
Features: Waterproof, Polycarbonate Case,
Stainless Steel Probe, On/Off Switch

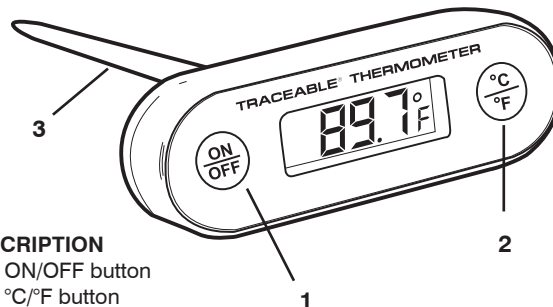

DESCRIPTION
1 - ON/OFF button
2 - °C/°F button
3 - Probe

ULTRA™ THERMOMETER ACCURACY

Ultra™ thermometers are tested at selected test points to be within tighter than normal tolerances to assist in providing improved accuracy. Other points will not necessarily fall within the same accuracy as those found at the selected test points, but will be within an accuracy of $\pm 1.0^{\circ}\text{C}$.

OPERATION

1. Press "ON/OFF" button to turn on unit.
2. Press the "°C/°F" button to select Fahrenheit or Celsius.

ALL OPERATIONAL DIFFICULTIES

If this unit does not function properly for any reason, replace the battery with a new high-quality battery (see "Battery Replacement" section). Low battery power can occasionally cause any number of "apparent" operational difficulties. Replacing the battery with a new fresh battery will solve most difficulties.

BATTERY REPLACEMENT

Erratic readings, a faint display, or no display are all indicators that the battery must be replaced. To replace the battery, loosen the two screws at the rear of the case. Remove the face of the thermometer taking care not to disconnect the internal wiring. Replace exhausted battery with a new 1.5V battery. Close the cover by tightening the screws. Replacement battery Cat. No. 1102.

NOTE: When replacing battery compartment cover, carefully ensure each cover is properly in place and screws tightened to maintain the unit's water-resistance. Incorrectly installed batteries may damage electronics.

WARRANTY, SERVICE, OR RECALIBRATION

For warranty, service, or recalibration, contact:

CONTROL COMPANY

4455 Rex Road
Friendswood, Texas 77546 USA
Ph. 281-482-1714 Fax 281-482-9448
E-mail sales@control3.com
www.control3.com

Control Company is ISO 9001 Quality-Certified by DNV and ISO 17025 accredited as a Calibration Laboratory by A2LA.

Cat. No. 4370 / 4369
Traceable® Food/Waterproof/Piercing/Drop-Proof Thermometer Instructions
92-4370-00
Rev. 3 083006

Traceable® is a registered trademark of Control Company