
Technical Services

Technical Services provides information on Berkshire products, quality programs and services.

Berkshire's demanding QA program has been developed to evaluate the suitability of Berkshire products to perform various functions. In addition to guaranteeing consistent product quality, the data generated by these tests is vital to Berkshire's quality assurance and product development processes.

Berkshire's quality system is registered to the ISO 9001 standard. If you would like more information Regarding Berkshire's QA Program or Test Methods please contact a sales representative or the Customer Service Department.

Berkshire Corporation
21 River Street
Great Barrington, MA 01230
800-242-7000
www.berkshire.com

BCR® and Choice® Cleanroom Socks

Washing

BCR® Cleanroom and Choice® Cleanroom socks can be re-laundered 2-3 times. It is recommended to wash on a gentle cycle with no bleach.

Note: Washer load size should be based on the washer manufacturer's guidelines, considering the density of the material and the ability of the socks to circulate in the drum during the wash cycle.

Drying

The dryer should be set at a low temperature setting.

Launderability

BCR® and Choice® Cleanroom Socks are re-launderable, however, the frequency of reuse is dependent upon the individual wearer and the duration of wear. Visual inspection of the socks is necessary after each laundering to determine their acceptability for continued use.

Autoclaving

The recommended autoclave cycle is 121°C for 30 minutes. This recommendation only covers the conditions the socks can withstand and does not ensure sterility is achieved.

It is the responsibility of the end user to validate the autoclave cycle to ensure the required results can be consistently achieved (considering all parameters, including but not limited to load size and load position).

Berkshire

