

Technical Data

Glucose Cysteine Agar Base w/ Thiamine

M433

Glucose Cysteine Agar Base w/ Thiamine with added blood or haemoglobin or hemin is used for cultivation and enumeration of *Francisella tularensis (Pasteurella tularensis)*.

Composition**

Ingredients	Gms / Litre
Meat peptone	3.000
Papaic digest of soyabean meal	10.000
Sodium chloride	5.000
Cysteine hydrochloride	1.000
Dextrose	25.000
Thiamine	0.0005
Agar	14.000
Final pH (at 25°C)	6.9±0.2

^{**}Formula adjusted, standardized to suit performance parameters

Directions

Suspend 58 grams in 1000 ml distilled water. Heat to boiling to dissolve the medium completely. Sterilize by autoclaving at 15 lbs pressure (121°C) for 15 minutes. Cool to 45-50°C and aseptically add sterile packed erythrocytes at a final concentration of 2% or 4-5% defibrinated sheep/rabbit blood. Mix well and pour into sterile Petri plates.

Principle And Interpretation

Francisella tularensis , a gram-negative aerobic bacillus, is the etiological agent of tularemia, which is primarily a disease of wild animals that is perpetuated in nature by ectoparasites, contaminated environment, cannibalism and acute or chronic carriers. Biting and blood sucking insects serve as vectors (1). Francisella (formerly known as Pasteurella) cannot be cultured on ordinary medium but require a complex medium containing blood or tissue extracts, thiamine and cysteine (2, 3). Glucose Cysteine Agar Base w/ Thiamine when supplemented with blood / haemoglobin is recommended for cultivation and enumeration of F.tularensis (Pasteurella tularensis)(4).

Meat peptone and papaic digest of soyabean meal provide essential growth nutrients. Dextrose serves as an easily metabolisable carbohydrate source while sodium chloride maintains the osmotic balance. Thiamine and cysteine hydrochloride serves as growth factor promoters required for culturing *Pasteurella*. Minute droplet like colonies develops in 48 hours.

Quality Control

Appearance

Cream to yellow homogeneous free flowing powder

Gelling

Firm, comparable with 1.4% Agar gel.

Colour and Clarity of prepared medium

.Basal medium: Amber coloured, clear to slightly opalescent gel forms. On addition of 4-5% sterile defibrinated sheep/rabbit blood: cherry red coloured opaque gel forms in Petri plates.

Reaction

Reaction of 5.8% w/v aqueous solution at 25°C. pH: 6.9±0.2

Growth

pН

6.70-7.10

Cultural Response

M433: Cultural characteristics observed with added 4-5% defibrinated sheep blood after an incubation at 35-37°C for 48-72 hours in presence of 10% CO₂

Organism

HiMedia Laboratories Technical Data

Francisella tularensis ATCC luxuriant 29684

Storage and Shelf Life

Store below 30°C in tightly closed container and prepared medium at 2-8°C. Use before expiry period on the label.

Reference

- 1.Murray P. R., Baron J. H., Pfaller M. A., Jorgensen J. H. and Yolken R. H., (Ed.), 2003, Manual of Clinical Microbiology, 8th Ed., American Society for Microbiology, Washington, D.C.
- 2. Collee J. G., Marmion B. P., Fraser A. G., and Simmons A., (Eds.), Mackie and McCartney Practical Medical Microbiology, 1996, 14th Edition, Churchill Livingstone, New York.
- 3.Manual of Diagnostic Tests and Vaccine for Terrestrial Animals, 2004, 5th Edi, OIE World Organization for Animal Health. 4.Atlas R. M., 2004, Handbook of Microbiological Media, Lawrence C. Parks, (Ed.), 3rd Edition, CRC Press, pg. no 717.

Revision: 1 / 2011

CE

Disclaimer :

User must ensure suitability of the product(s) in their application prior to use. Products conform solely to the information contained in this and other related HiMedia™ publications. The information contained in this publication is based on our research and development work and is to the best of our knowledge true and accurate. HiMedia™ Laboratories Pvt Ltd reserves the right to make changes to specifications and information related to the products at any time. Products are not intended for human or animal or therapeutic use but for laboratory, diagnostic, research or further manufacturing use only, unless otherwise specified. Statements contained herein should not be considered as a warranty of any kind, expressed or implied, and no liability is accepted for infringement of any patents.