


HOW TO USE THE LiveSlide®

Using the LiveSlide® microscope slide is easy. Simply follow these directions.

1. A LiveSlide® set contains 3 pieces, the base slide and a set of coverslips, all of which are reusable.
2. The 2 coverslips are different thicknesses. One is very thin, perfect for high magnification. The second is a bit thicker for all applications.
3. The coverslips have a thin, plastic film on both sides for protection – clear and green. Remove the films by peeling the film off or use a small piece of tape on the corner and pull back the film. Discard the film.

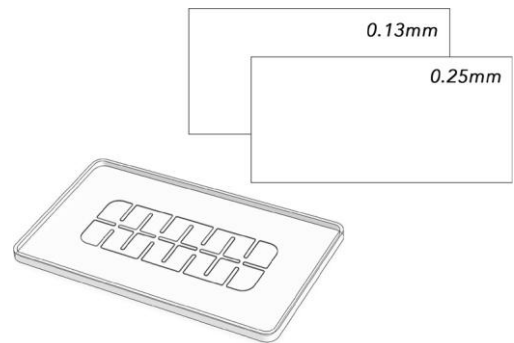

4. Gather your collected specimens.



5. Next, get the base slide and place 6-10 drops on it.



6. Gently place the coverslip over the culture well.



7. Place the LiveSlide® on a microscope stage or under a digital microscope and enjoy exploring the magnified world.

8. Rinse with warm water, and, when dry, store for future use.