

# ABSOLUTE ZERO® DRY ICE SNOW MAKER

## Instruction Manual




Innovative shipping and  
packaging solutions by

***Polar Tech***  
INDUSTRIES, INC.


# Absolute Zero® Dry Ice Snow Maker

## Installation Instructions


- 1 Connect female end of hose to snow maker.  
(Ensure connection is tight)

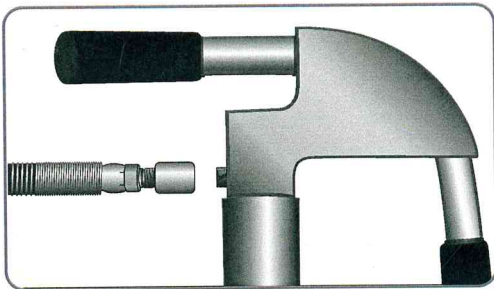


- 2 Connect male end of hose to brass tee.  
(Ensure connection is tight)

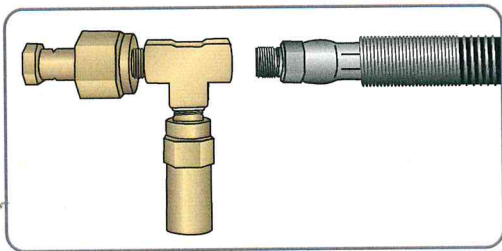


- 3 Insert washer into tank connector.

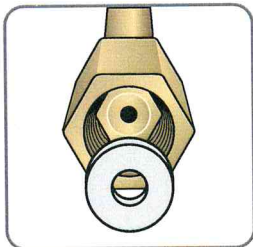


- 4** Connect tank connector to “liquid” outlet on CO<sub>2</sub> tank.

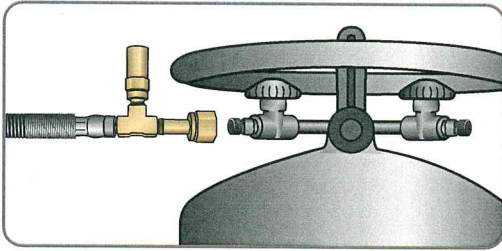

- 5** Open tank valve 1/2 to 3/4 of a turn

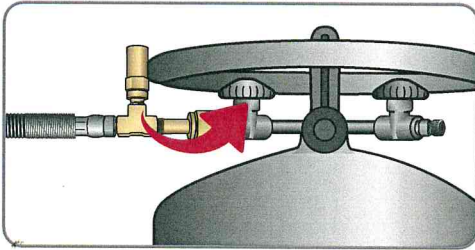

- 6** Fill container with snow.

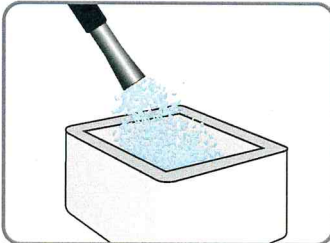



# WARNING

- Dry Ice is extremely cold ( $-109.3^{\circ}\text{F}$ ,  $-78.5^{\circ}\text{C}$ ) and, if not properly handled, can cause severe cold burns and frostbite.
- Do not touch Dry Ice with bare hands, use heavy insulated gloves. Rubber and plastic gloves do not provide adequate protection.
- Use protective eyewear when making or handling Dry Ice.
- Dry Ice is not edible. Do not put Dry Ice in your mouth and do not allow anyone else to do so.
- Keep Dry Ice out of the reach of children and animals.
- Do not put Dry Ice in any tightly closed container.
- Do not add Dry Ice to any liquid or any liquid to Dry Ice.
- Make and use Dry Ice only in a well-ventilated area.
- Before operating the Dry Ice Snow Maker, make sure all tank and hose connections are tight. All connections must be tightened with a wrench.
- Consult your  $\text{CO}_2$  tank supplier about recommended procedures for proper tank installation, storage and operation.

**Polar Tech Industries, Inc.**

1-800-ICE-BRIX

[www.polar-tech.com](http://www.polar-tech.com)

Copyright © 2012 Polar Tech Industries, Inc.