

Bransonic® Ultrasonic Cleaning Bath Accessories



General Description

Bransonic accessories are designed to maximize the effectiveness of your Bransonic baths. It is important that parts do not rest on the bottom of the bath, as it can inhibit the ultrasonic activity and possibly damage the unit.

A variety of perforated trays, mesh baskets, beakers, and support racks give you the option of direct or indirect cleaning. All five models come with covers to allow the cleaning bath to heat up more quickly.

Direct Cleaning Method

- Suspends parts using a perforated tray, mesh basket, or a support rack.
- Permits sound waves to penetrate directly to the parts.

Indirect Cleaning Method

- Suspends parts using a solid tray, or beakers with a beaker cover.
- Ideal for cleaning small parts.
- Discarded soil cleans easily from a beaker or solid tray.
- Indirect method is ideal for sample preparation.






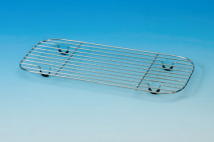

Consult with your Branson representative about the Bransonic accessories and solution best suited for your application needs. **Visit us at www.Bransonic.com**

Additional Features

- Perforated trays, solid insert trays, and wire mesh baskets have built-in handles specially designed to prevent accessories from touching the tank bottom.
- Support racks allow total volume of the tank to be used while keeping parts, fixtures, or containers about ½" above the tank bottom.
- Support racks also are useful as countertop drain racks, holding parts above the countertop surface.

BRANSON

Bransonic Accessories – Order using part numbers listed under model numbers

	Ordering Part No.	Model 1800	Model 2800	Model 3800	Model 5800	Model 8800
 Cover (Supplied with Unit)		100-032-515	100-032-517	100-032-519	100-032-510	000-917-050 (SS)
 *Perforated Tray		100-410-160	100-410-162	100-410-164	100-410-166	100-410-168
 Mesh Basket		100-916-333	100-916-334	100-916-335	—	—
 Solid Insert Tray		100-410-170	100-410-172	100-410-174	100-410-176	100-410-178
 Beaker Positioning Cover		(1x600 ml) 100-032-516	(2x600 ml) 100-032-518 (2x250 ml) 100-032-527	(2x600 ml) 100-032-520 (3x250 ml) 100-032-528	(4x600 ml) 100-032-511	(6x600 ml) CPN-246-014 (SS)
 Support Racks		CPN-916-039	CPN-916-040	CPN-916-041	CPN-916-042	CPN-916-043
 Beakers (All Models)		250 ml Glass – 000-140-001 400 ml PP – 000-265-061 600 ml Glass – 000-140-004 600 ml SS – 000-410-055				

*Note: Models 1800, 2800 and 3800: Perforated holes run lengthwise towards the bottom; 3 rows only.

Americas

Branson Ultrasonics Corp.
41 Eagle Road
Danbury, CT 06810, USA
T: 203-796-0400
F: 203-796-2250
Toll Free: 800-732-9262
www.bransonultrasonics.com

Europe and Africa and Middle East

Branson Ultrasonic SA
Ch. du Faubourg-de-Cruiseilles 9
1227 Carouge
Switzerland
T: +41(0)22 3048353
F: +41(0)22 3048359
www.branson.ch

Asia

Branson Ultrasonics (Shanghai) Co., Ltd.
758 Rong Le Dong Road
Song Jiang, Shanghai, PRC, 201613
T: 86-21-3781-0588
F: 86-21-5774-5100
www.branson.com.cn

All specifications subject to change without notice. All dimensions are nominal.



© Branson Ultrasonics Corporation 2016.
The Emerson logo is a trademark and service mark
of Emerson Electric Co.
Printed in the U.S.A.

BCL-0004-12